



**DAICEL CHIRAL TECHNOLOGIES
(INDIA) PRIVATE LIMITED**

SUBSIDIARY OF DAICEL CHEMICAL INDUSTRIES, LTD.

IPC-USP 8th ASM 2009

“Chiral Impurity Methods – Case Study”

**Ch. LAKSHMI NARAYANA
DAICEL CHIRAL TECHNOLOGIES (INDIA) PVT LTD**





DAICEL CHIRAL TECHNOLOGIES
(INDIA) PRIVATE LIMITED

SUBSIDIARY OF DAICEL CHEMICAL INDUSTRIES, LTD.

Overview

- Introduction: Chiral Impurity methods
- Types of Chiral Stationary Phases (CSPs) for HPLC
 - Polysaccharide-CSPs
 - Coated CSPs
 - immobilised CSPs
- Case study:
 - Oxaliplatin Chiral HPLC method
 - Lamivudine Chiral HPLC method
 - Clopidogrel Bisulphate Chiral HPLC method





DAICEL CHIRAL TECHNOLOGIES
(INDIA) PRIVATE LIMITED

SUBSIDIARY OF  DAICEL CHEMICAL INDUSTRIES, LTD.

Chiral Impurity Methods – Introduction

- Chiral GC
- Chiral HPLC
- Chiral HPCE
- Chiral SFC





DAICEL CHIRAL TECHNOLOGIES
(INDIA) PRIVATE LIMITED

SUBSIDIARY OF DAICEL CHEMICAL INDUSTRIES, LTD.

Chiral Impurity Methods – Types of CSP

Types of CSPs and their loading capacities

Type	CSPs	Loading capacity (mg solute / g CSP)
I	Pirkle type (Brush type)	1-50
II	Polysaccharide derivatives	5-150
III	Macrocyclic type	
	Cyclodextrins	0.1-5
	Glycopeptides	0.1-5
	Chiral Crown ether	0.1-5
IV	Ligand exchange	0.1-1
V	Protein type	0.1-0.2





DAICEL CHIRAL TECHNOLOGIES
(INDIA) PRIVATE LIMITED

SUBSIDIARY OF DAICEL CHEMICAL INDUSTRIES, LTD.

Chiral Impurity Methods – Types of CSP


Which type of CSP ?

1999

2001

2003

2005

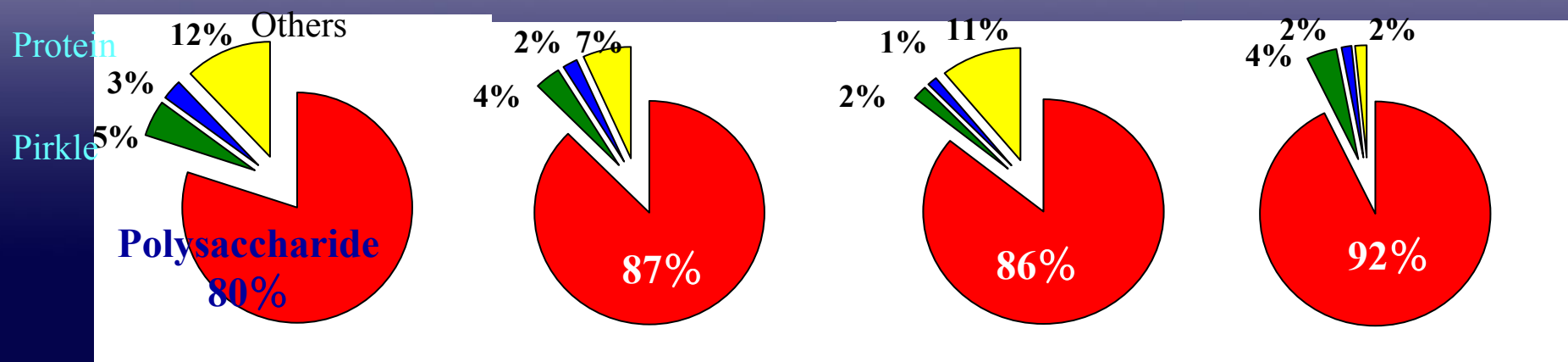

1995-2003 : *Tetrahedron Asymmetry*

2005 : *J. Am. Chem. Soc.*






DAICEL CHIRAL TECHNOLOGIES
(INDIA) PRIVATE LIMITED

SUBSIDIARY OF DAICEL CHEMICAL INDUSTRIES, LTD.


Chiral Impurity Methods – Types of CSP

Coated polysaccharide-derived CSPs

Cellulose derivatives

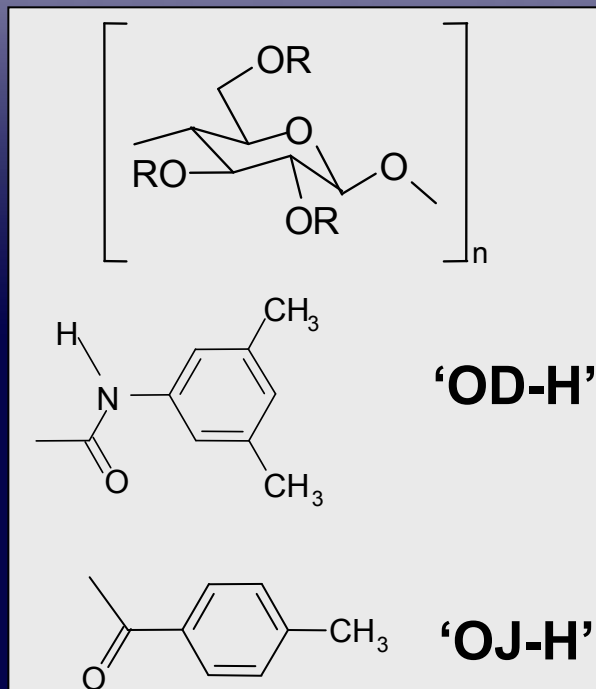

Amylose derivatives

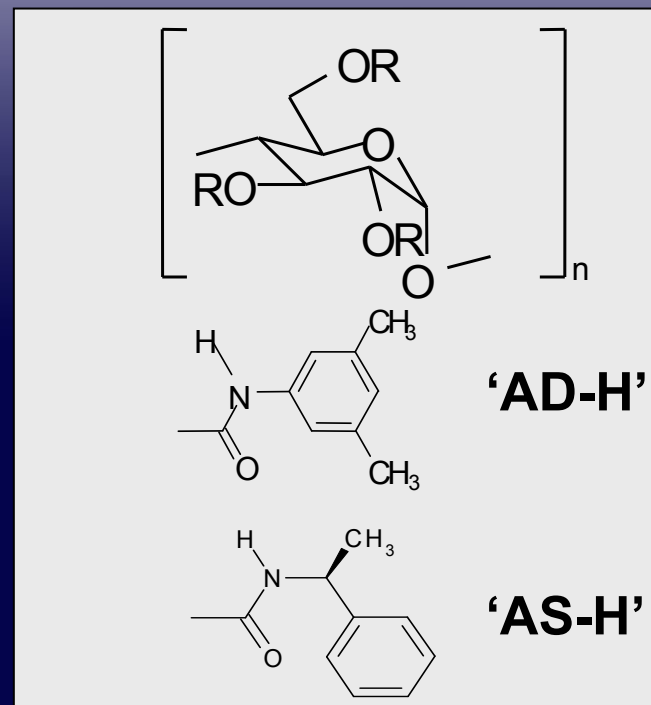




DAICEL CHIRAL TECHNOLOGIES
(INDIA) PRIVATE LIMITED


SUBSIDIARY OF DAICEL CHEMICAL INDUSTRIES, LTD.

Chiral Impurity Methods – Types of CSPs


Immobilised polysaccharide-derived CSPs

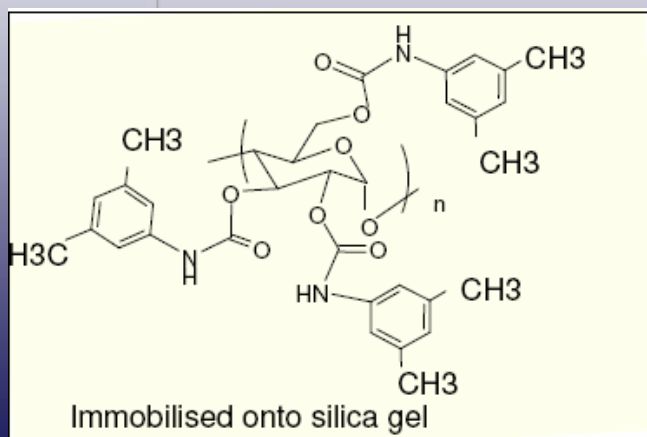

CHIRALPAK IA

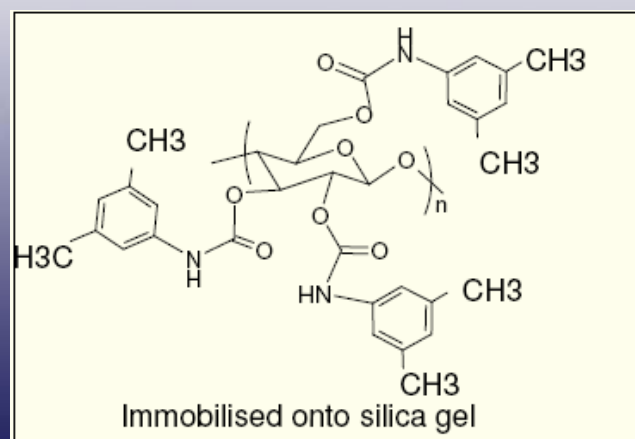

CHIRALPAK IB

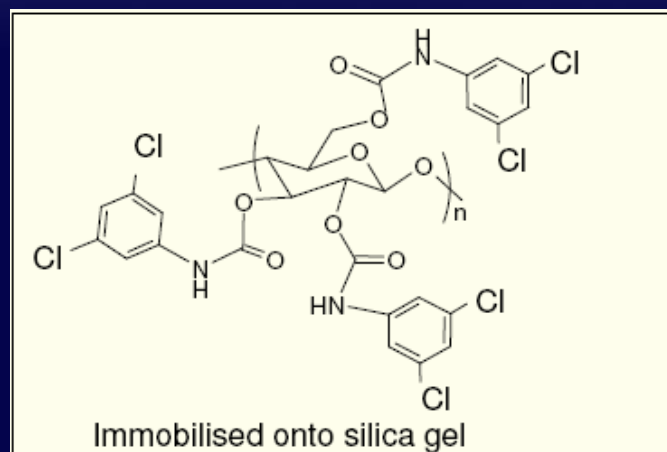

CHIRALPAK IC







DAICEL CHIRAL TECHNOLOGIES
(INDIA) PRIVATE LIMITED

SUBSIDIARY OF DAICEL CHEMICAL INDUSTRIES, LTD.

Chiral Impurity Methods-Case Study

- Case study-1:
Oxaliplatin, Chemotherapy drug
- Case study-2:
Lamivudine, Anti retroviral drug
- Case study-3:
Clopidogrel Bisulphate, Anti platelet drug

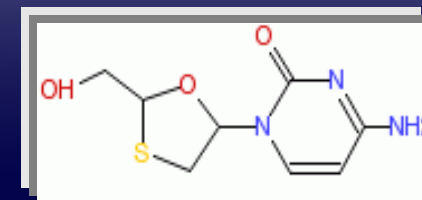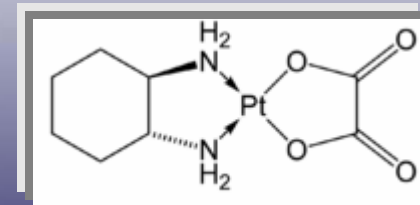




DAICEL CHIRAL TECHNOLOGIES
(INDIA) PRIVATE LIMITED

SUBSIDIARY OF DAICEL CHEMICAL INDUSTRIES, LTD.

Chiral Impurity Methods – Case Study 1

Oxaliplatin

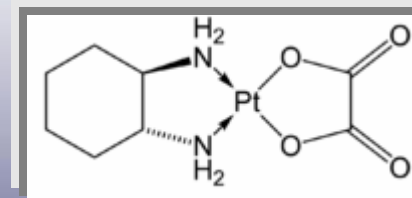

Pharmacopeial method conditions:

Column: CHIRALCEL OC 4.6 x 250mm

Mobile Phase: Ethanol / MeOH (30/70)


Flow Rate: 0.3 mL/min; Injection Volume: 20 μ L

Detection : 254 nm by UV; Temp: 40° C

Concentration: 0.6 mg/mL

R_s is NLT 1.5

Limit: 0.1%

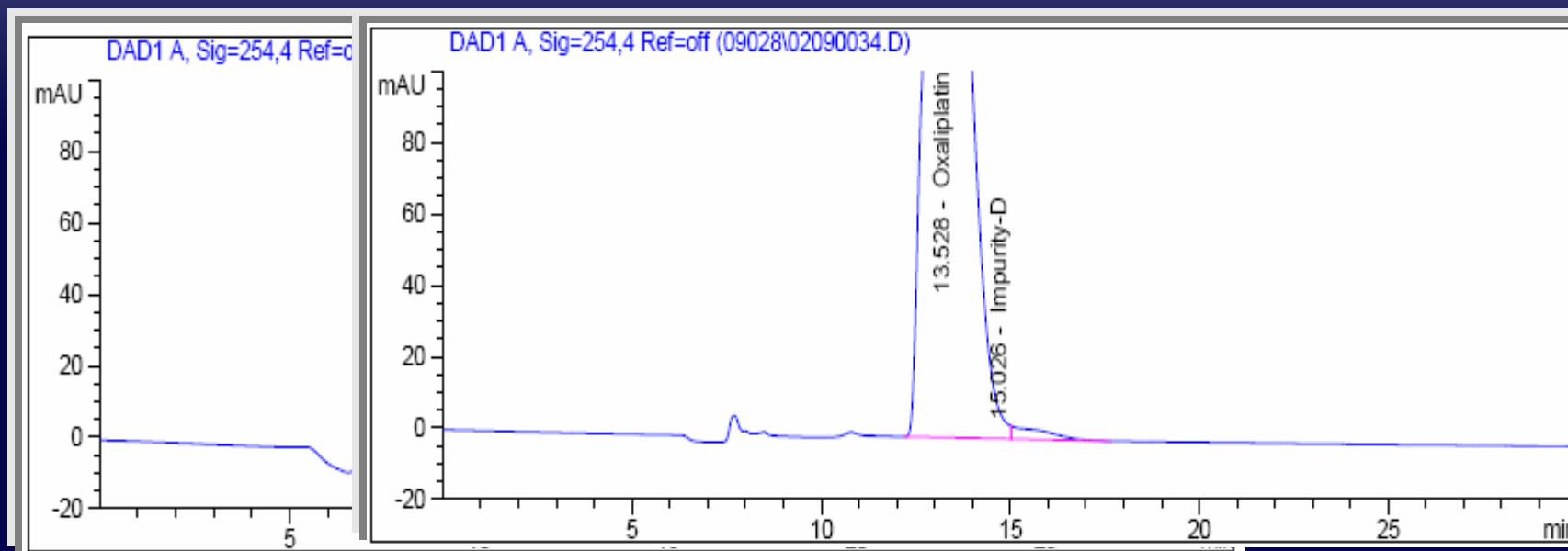



DAICEL CHIRAL TECHNOLOGIES
(INDIA) PRIVATE LIMITED

SUBSIDIARY OF DAICEL CHEMICAL INDUSTRIES, LTD.

Chiral Impurity Methods – Case Study 1

Oxaliplatin

Cost-effective method conditions:

Column: CHIRALPAK IC 4.6 x 250mm

Mobile Phase: Ethanol / MeOH (30/70)


Flow Rate: 1.0 mL/min; Injection Volume: 20 μ L

Detection : 254 nm by UV; Temp: 40° C

Concentration: 0.6 mg/mL

$R_s > 9.0$

LOQ: 0.03%

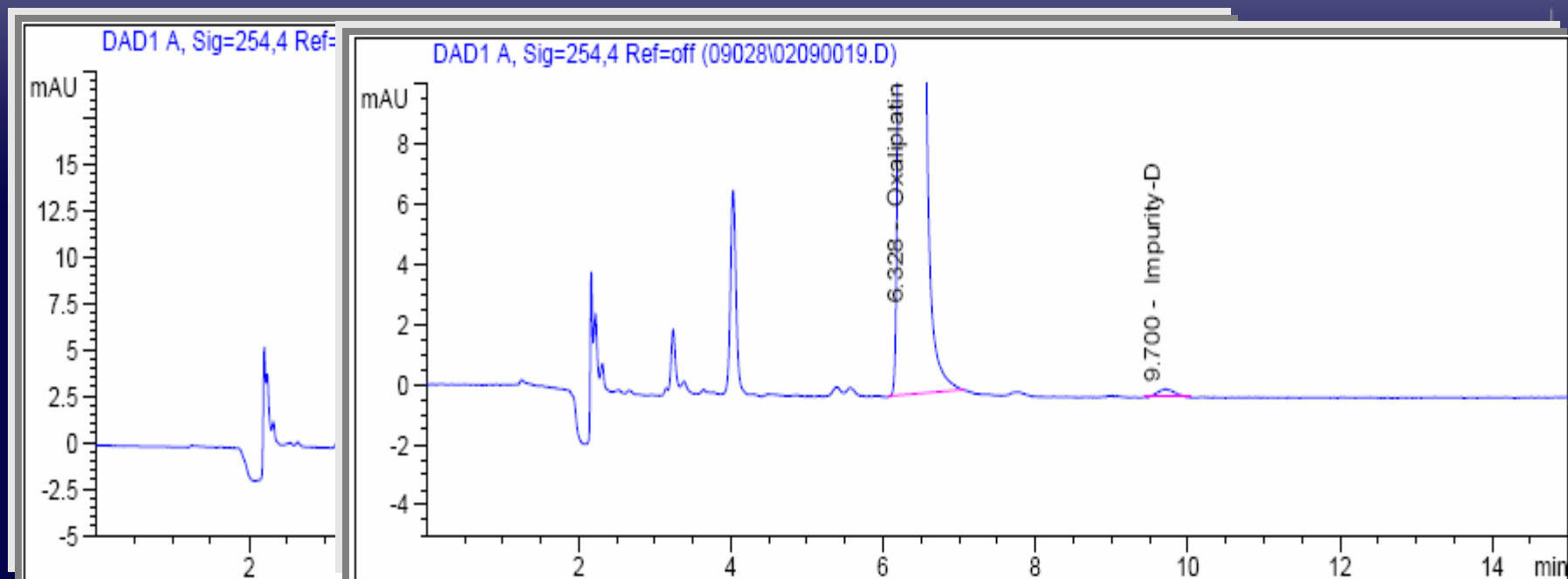



DAICEL CHIRAL TECHNOLOGIES
(INDIA) PRIVATE LIMITED

SUBSIDIARY OF DAICEL CHEMICAL INDUSTRIES, LTD.

Chiral Impurity Methods – Case Study 1

Oxaliplatin

Cost-effective method conditions:


Column: CHIRALPAK IC 4.6 x 250mm

Mobile Phase: Ethanol / MeOH (30/70)

Flow Rate: 1.0 mL/min; Injection Volume: 20 μ L

Detection : 254 nm by UV; Temp: 40° C

Concentration: 0.6 mg/mL

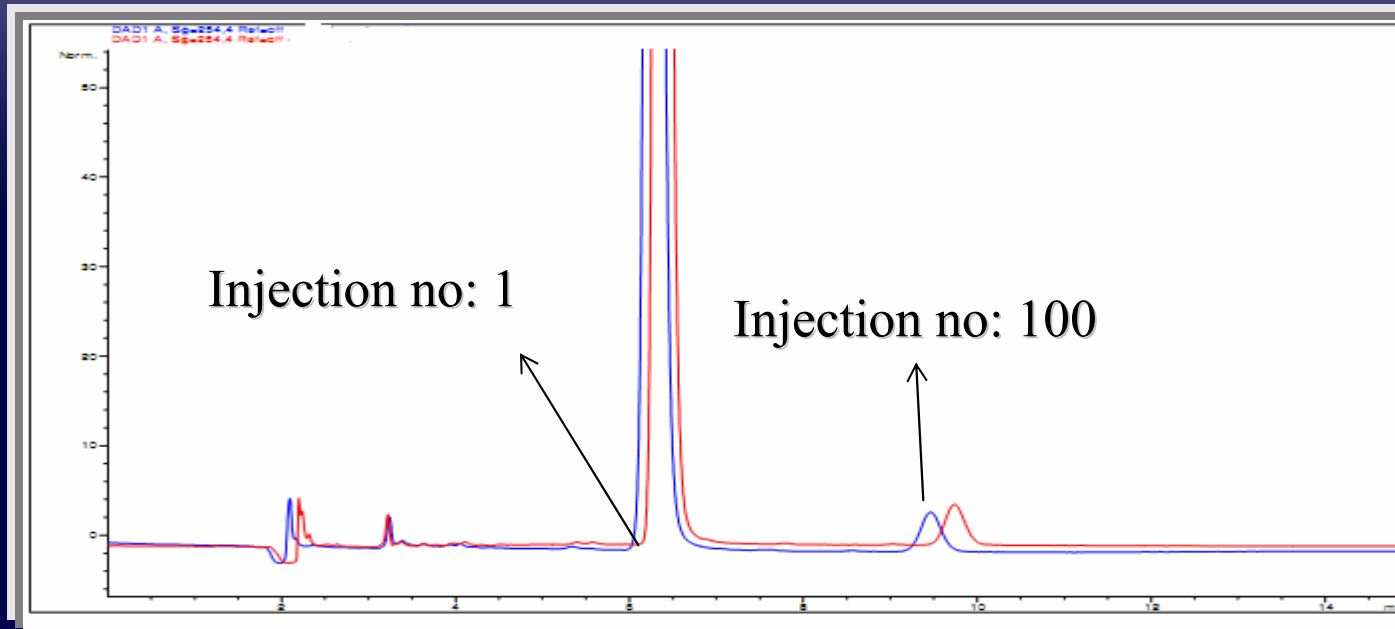




DAICEL CHIRAL TECHNOLOGIES
(INDIA) PRIVATE LIMITED

SUBSIDIARY OF DAICEL CHEMICAL INDUSTRIES, LTD.

Chiral Impurity Methods – Case Study 2

Lamivudine

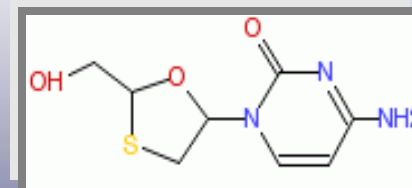

Pharmacopeial method conditions:

Column: L45 (Cyclobond I 2000 SP, 4.6 x 250mm, 5 μ)


Mobile Phase: 0.1M Ammonium acetate/MeOH (95/05)

Flow Rate: 1.0 mL/min; **Conc:** 0.25 mg / mL; **Inj Vol:** 10 μ L

Detection : 270 nm by UV; **Temperature:** 25 ° C

R_s is NLT 1.5

Impurity limit: 0.3%

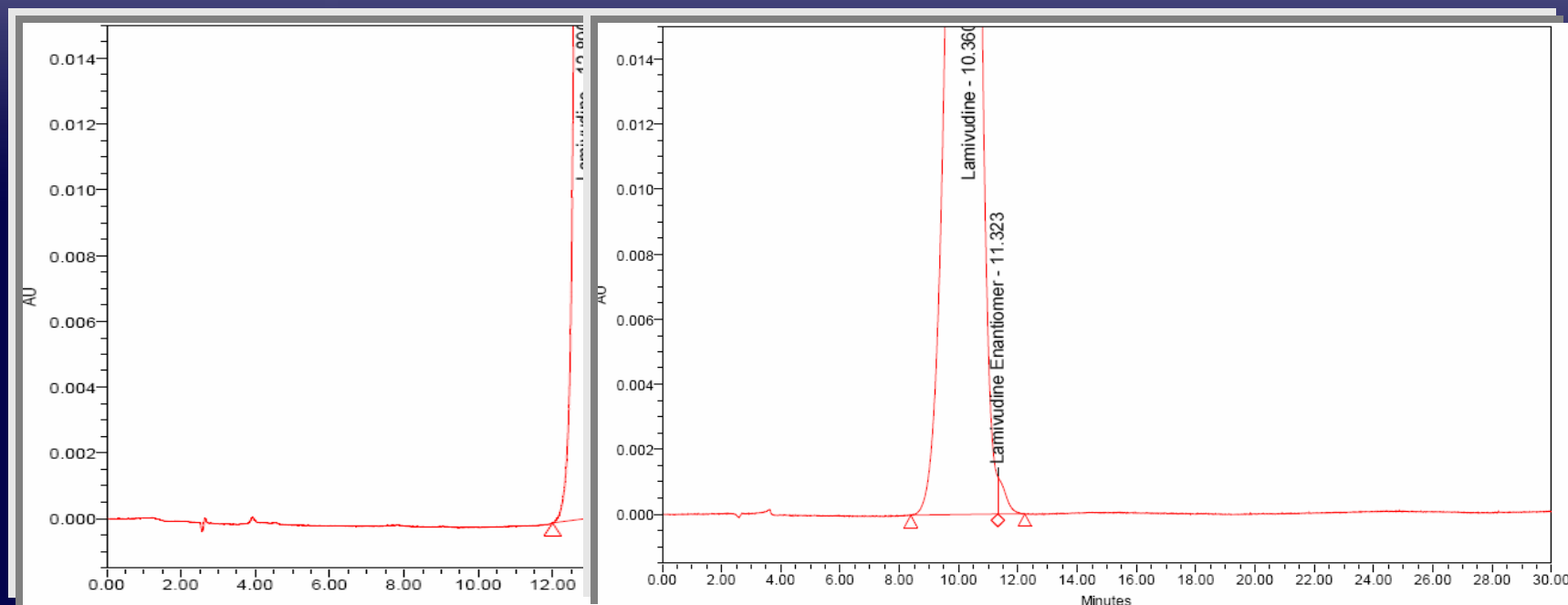



DAICEL CHIRAL TECHNOLOGIES
(INDIA) PRIVATE LIMITED
SUBSIDIARY OF DAICEL CHEMICAL INDUSTRIES, LTD.

Chiral Impurity Methods – Case Study 2

Lamivudine


Cost-effective method conditions:

Column: CHIRALPAK IC (4.6 x 250) mm, 5 micron

Mobile Phase: Ethanol/2-Propanol/DEA (90/10/0.1, v/v/v)

Flow Rate: 0.5 mL/min; Conc: 0.25 mg/mL; Inj Vol: 10 μ L

Detection : 270 nm by UV; Temperature: 25 ° C; Diluent : MP

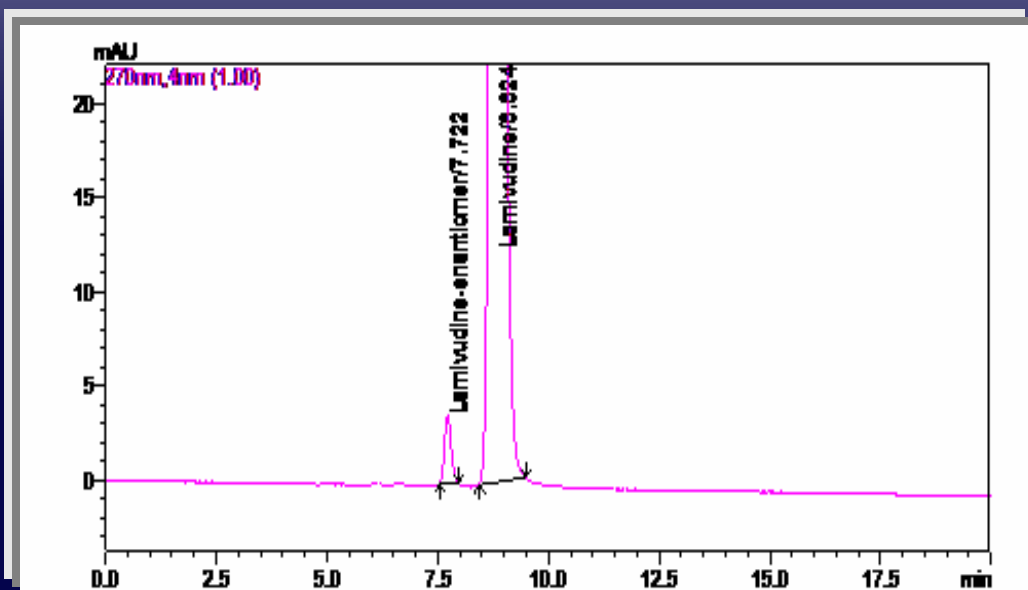

$R_s > 3.0$

LOQ: 0.1%





DAICEL CHIRAL TECHNOLOGIES
(INDIA) PRIVATE LIMITED
SUBSIDIARY OF DAICEL CHEMICAL INDUSTRIES, LTD.

Chiral Impurity Methods – Case Study 2

Lamivudine


Cost-effective method conditions:

Column: CHIRALPAK IC (4.6 x 250) mm, 5 micron

Mobile Phase: Ethanol/2-Propanol/DEA (90/10/0.1, v/v/v)

Flow Rate: 0.5 mL/min

Detection : 270 nm by UV; Temperature: 25 ° C

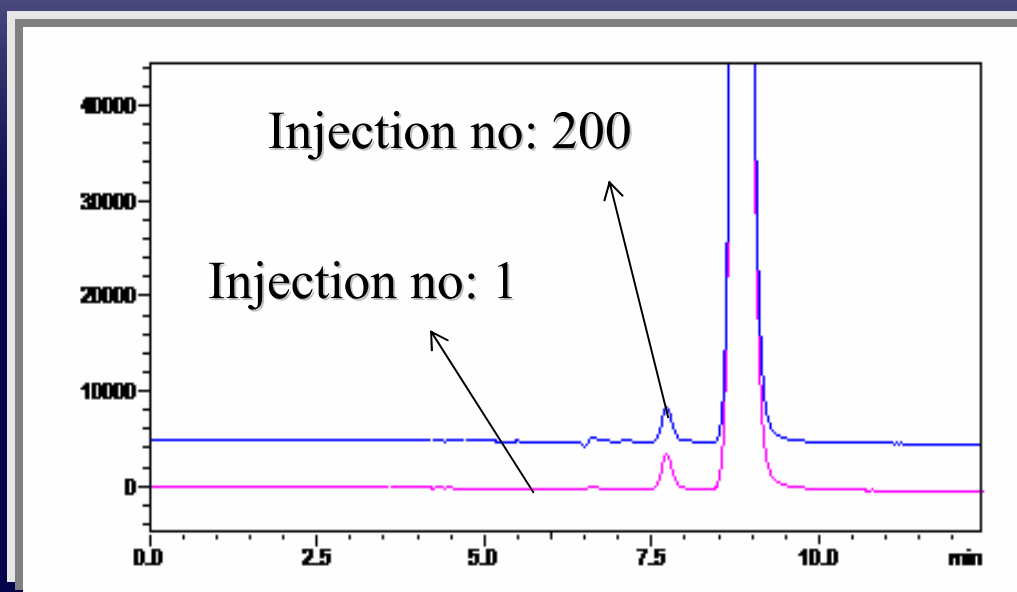



DAICEL CHIRAL TECHNOLOGIES
(INDIA) PRIVATE LIMITED

SUBSIDIARY OF DAICEL CHEMICAL INDUSTRIES, LTD.

Chiral Impurity Methods – Case Study 3

Clopidogrel bisulphate


Pharmacopeial method conditions:

Column: L 57 (Ultron ES-OVM (4.6 x 150) mm)

Mobile Phase: 10mM Phosphate buffer/ACN (75/25)

Flow Rate: 1.0 mL/min; Conc: 0.5 mg/mL

Detection : 220 nm by UV

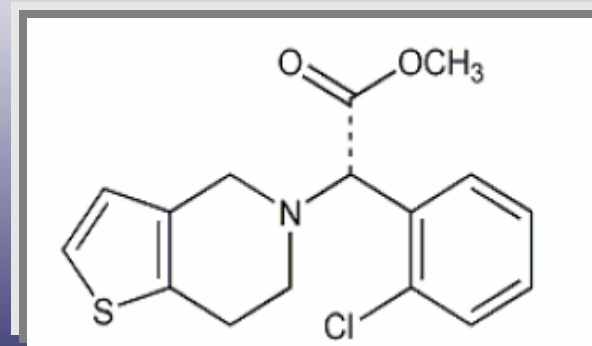

- Method objective:
 1. Clopidogrel bisulphate related substances quantification including chiral impurity
 2. Clopidogrel bisulphate assay determination






DAICEL CHIRAL TECHNOLOGIES
(INDIA) PRIVATE LIMITED


SUBSIDIARY OF DAICEL CHEMICAL INDUSTRIES, LTD.

Chiral Impurity Methods – Case Study 3


Clopidogrel and its related compounds

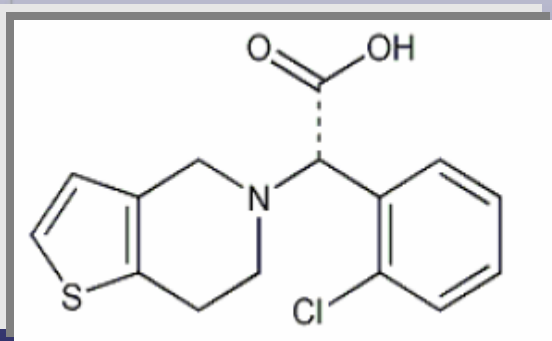

Imp-A
Limit: 0.2%

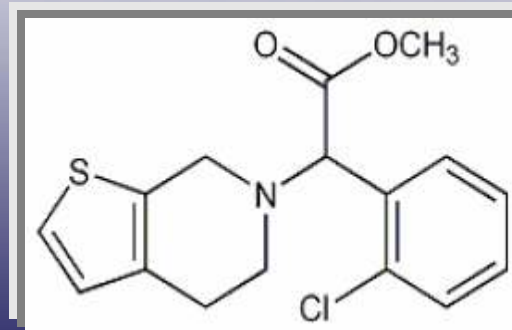

Imp-B
Limit: 0.3%

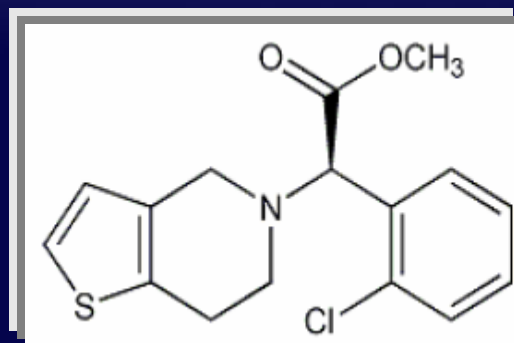

Imp-C
Limit: 1.0%





DAICEL CHIRAL TECHNOLOGIES
(INDIA) PRIVATE LIMITED


SUBSIDIARY OF DAICEL CHEMICAL INDUSTRIES, LTD.

Chiral Impurity Methods – Case Study 3

Clopidogrel bisulphate

Pharmacopeial method

SS criteria: Rs between Imp B1 and Clopidogrel NLT 2.5

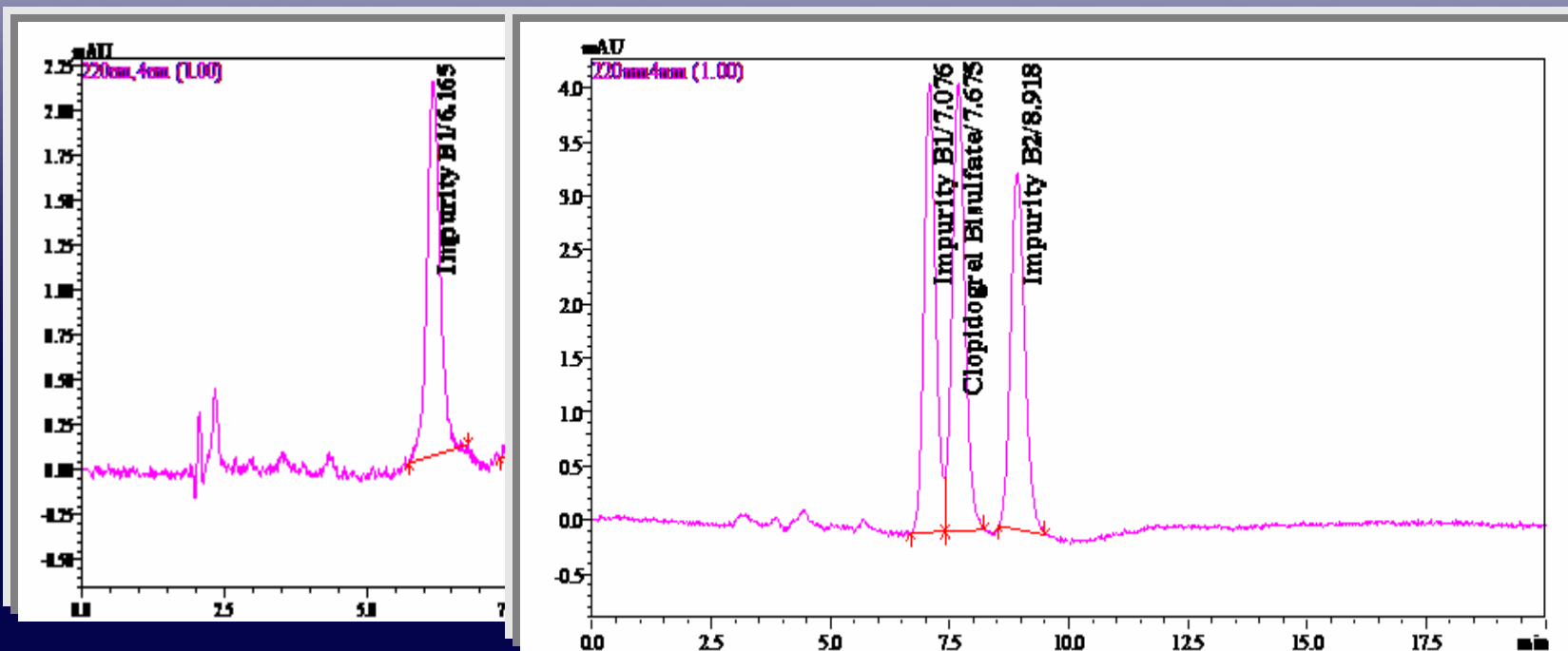




DAICEL CHIRAL TECHNOLOGIES
(INDIA) PRIVATE LIMITED

SUBSIDIARY OF DAICEL CHEMICAL INDUSTRIES, LTD.

Chiral Impurity Methods – Case Study 3

Clopidogrel bisulphate

Pharmacopeial method

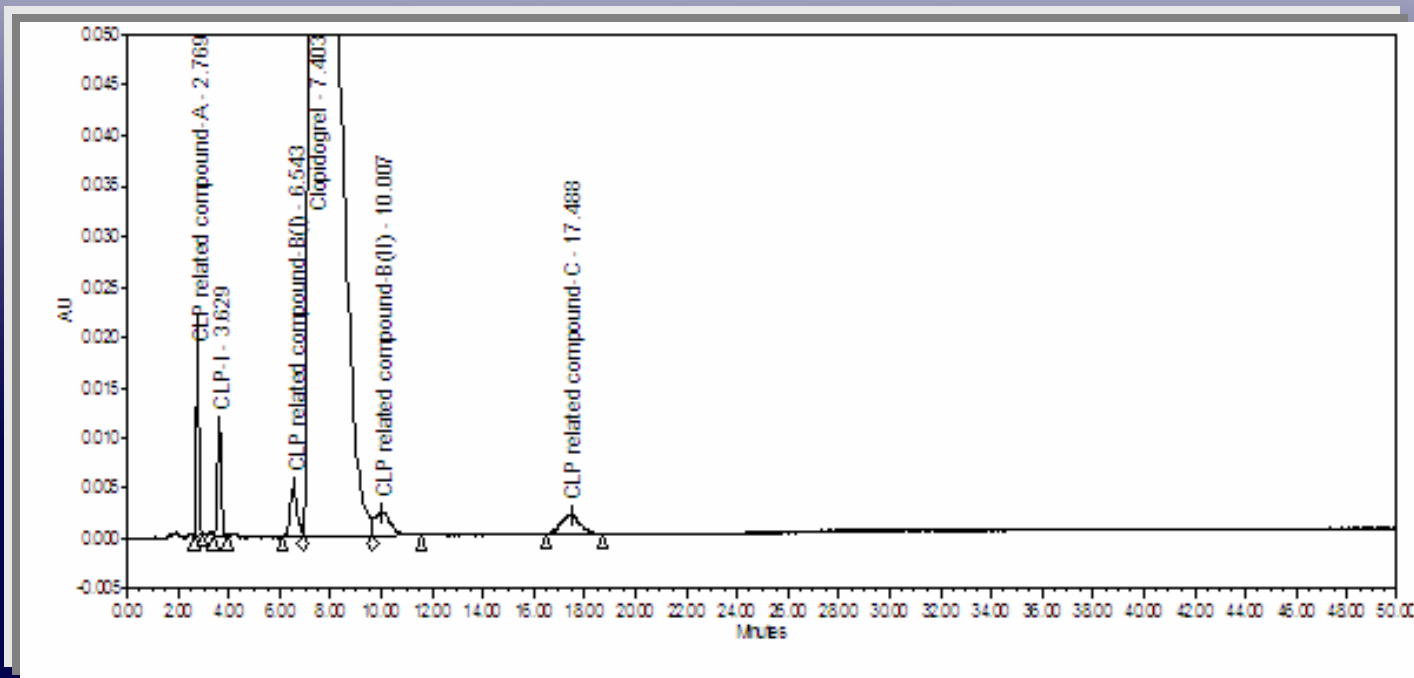



DAICEL CHIRAL TECHNOLOGIES
(INDIA) PRIVATE LIMITED

SUBSIDIARY OF DAICEL CHEMICAL INDUSTRIES, LTD.

Chiral Impurity Methods – Case Study 3

Clopidogrel bisulphate


Cost effective method conditions:

Column: CHIRALCEL OJ-H (4.6 x 250) mm, 5 micron

Mobile Phase: Methanol/DEA (100/0.1, v/v)

Flow Rate: 1.0 mL/min; Conc: 0.5 mg/mL; Diluent : Ethanol

Detection : 220 nm by UV; Temperature: 25 ° C

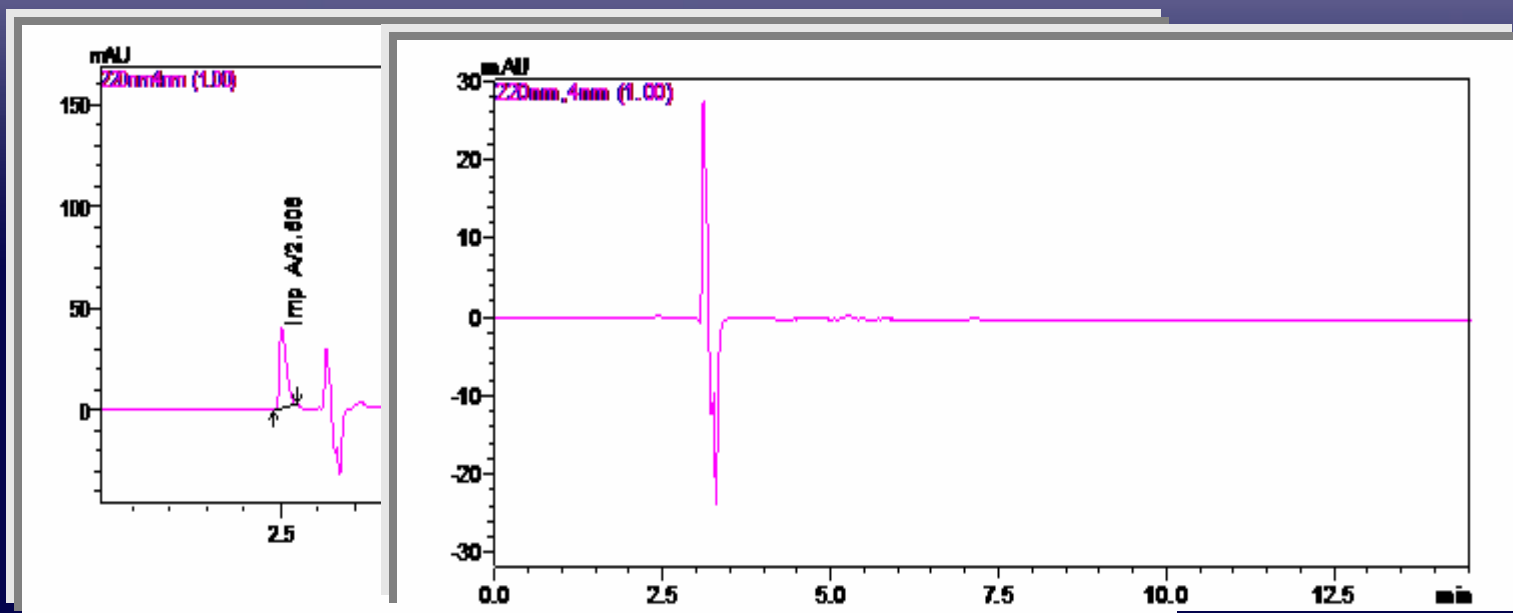



DAICEL CHIRAL TECHNOLOGIES
(INDIA) PRIVATE LIMITED

SUBSIDIARY OF DAICEL CHEMICAL INDUSTRIES, LTD.

Chiral Impurity Methods – Case Study 3

Clopidogrel bisulphate


Cost-effective method conditions:

Column: CHIRALCEL OJ-H (4.6 x 250) mm, 5 micron

Mobile Phase: Methanol/DEA (100/0.1, v/v)

Flow Rate: 1.0 mL/min; Diluent : Ethanol

Detection : 220 nm by UV; Temperature: 25 ° C

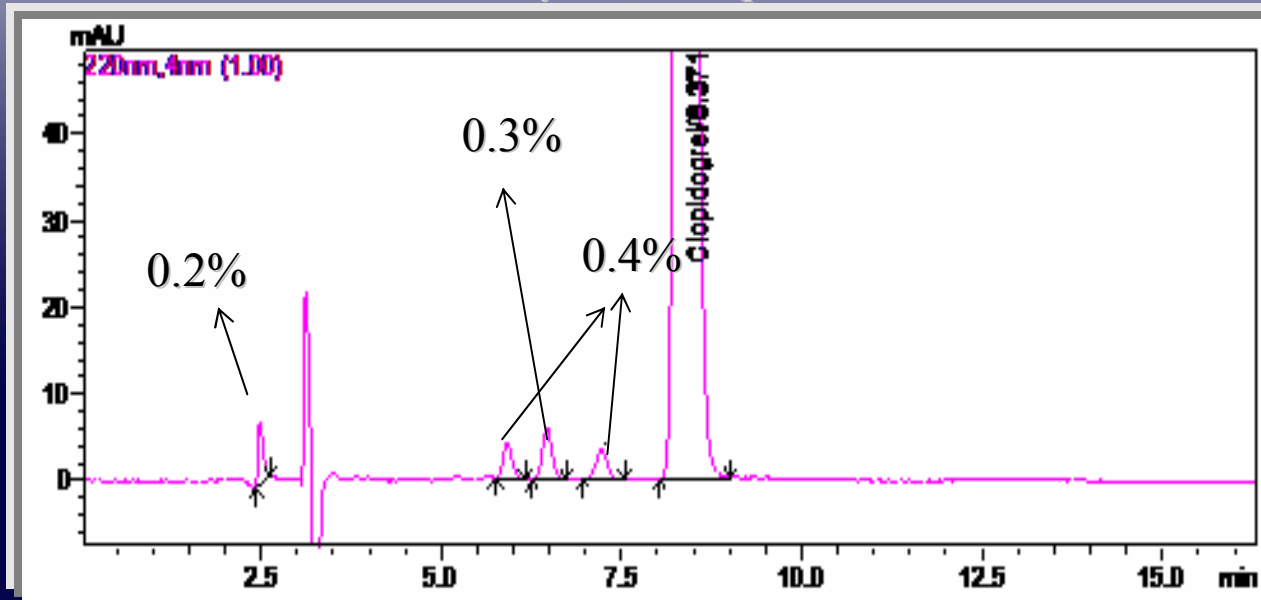



DAICEL CHIRAL TECHNOLOGIES
(INDIA) PRIVATE LIMITED

SUBSIDIARY OF DAICEL CHEMICAL INDUSTRIES, LTD.

Chiral Impurity Methods – Case Study 3

Clopidogrel bisulphate


Cost-effective method conditions:

Column: CHIRALCEL OJ-H (4.6 x 250) mm, 5 micron

Mobile Phase: Methanol/DEA (100/0.1, v/v)

Flow Rate: 1.0 mL/min; Diluent : Ethanol

Detection : 220 nm by UV; Temperature: 25 ° C

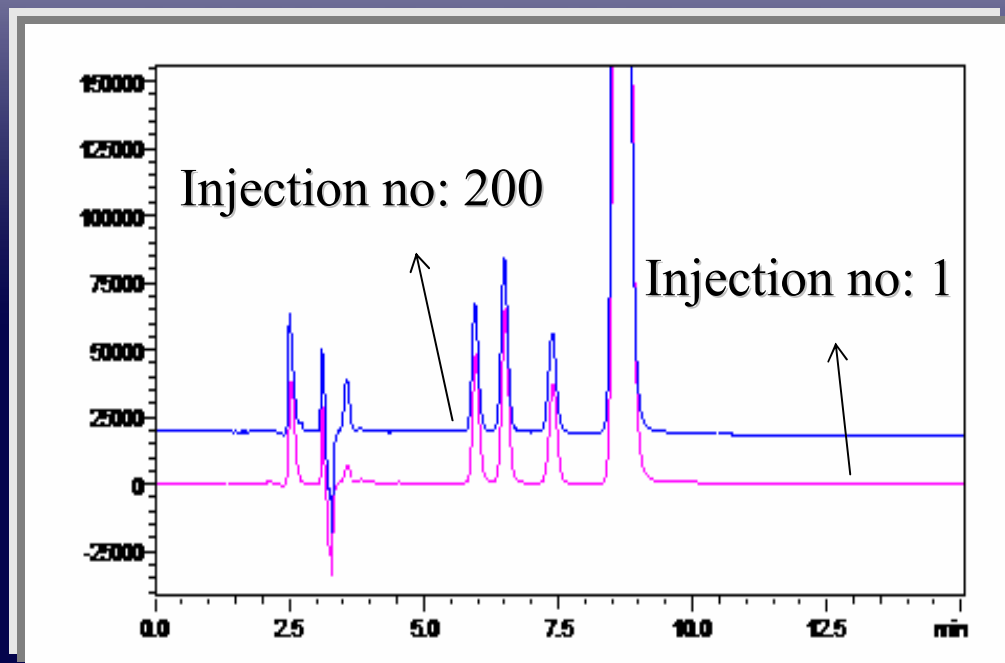



DAICEL CHIRAL TECHNOLOGIES
(INDIA) PRIVATE LIMITED
SUBSIDIARY OF DAICEL CHEMICAL INDUSTRIES, LTD.

Chiral Impurity Methods – Case Study 3

Clopidogrel bisulphate

Chromatographic Characteristics

Compound peak	Rs	N	T
Imp A	-	3907	1.4
Imp B1	17.4	11520	1.3
Imp C	2.3	11554	1.1
Imp B2	2.9	11063	1.1
Clopidogrel	3.6	10235	1.3

Note: Column memory effects may influence the elution of Imp A, hence recommended to dedicate the column for this application.





DAICEL CHIRAL TECHNOLOGIES
(INDIA) PRIVATE LIMITED

SUBSIDIARY OF DAICEL CHEMICAL INDUSTRIES, LTD.

Conclusions

- Chiral HPLC using polysaccharide derived CSPs is a versatile tool to estimate chiral impurities in drug substances .
- Perhaps, the chiral impurity methods described in Pharmacopeia for Oxaliplatin, Lamivudine & Clopidogrel bisulphate are utilising the columns, which are not so durable and hence results into high analytical cost.
- It would be beneficial to adapt new generation chiral column chemistry available today to develop an efficient & cost effective chiral impurity methods.





DAICEL CHIRAL TECHNOLOGIES
(INDIA) PRIVATE LIMITED

SUBSIDIARY OF DAICEL CHEMICAL INDUSTRIES, LTD.

Acknowledgments

- Mr. Raghuram, General Manager, Hetero Drugs Ltd for his continuous co-operation during Lamivudine method development
- My Colleagues Mr. Thirupathi & Mr. Srinivasu for their significant contribution during method development
- USP India for the lecture opportunity





DAICEL CHIRAL TECHNOLOGIES
(INDIA) PRIVATE LIMITED

SUBSIDIARY OF  **DAICEL CHEMICAL INDUSTRIES, LTD.**

THANK YOU ALL

